

## Fractional Efficiency Measurement System

## FMS 375



Fractional efficiency measuring system FMS 375.

The fractional efficiency measurement system FMS 375 bundles core competences of Topas in one compact and robust instrument for highly-resolved aerosol analyses regarding particle size and concentration.

The FMS 375 enables an automated aerosol sampling as well as excess air refeeding via a switching unit, aerosol conditioning by two static dilution systems and a precise aerosol characterisation based on the aerosol spectrometer LAP 325.

### Applications


- characterisation of particle size distributions and particle numbers of aerosols
- evaluation of fractional separation efficiency of oil mist separators and filters
- compact design for use at engine test stations or in laboratory test setups (e.g. SPT 140)

### Features

- operable at both over and lower pressure
- automated and remote-controlled operation
- mobile use (optional trolley kit)

### Principle of operation

The FMS 375 has to be connected to the sampling points of a test set-up. These are then individually controlled via the automated switching unit. If necessary, the sample is conditioned before characterisation by means of the dilution systems. The dilution factor can be varied in three steps. The particle inlet concentration defines the necessity for none, one or two dilution systems for lowering the concentration towards the acceptable concentration range for the aerosol spectrometer.

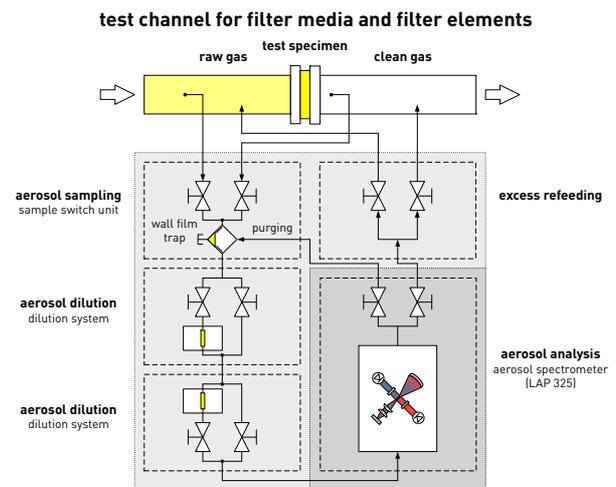

Principle of operation of the fractional efficiency measurement system FMS 375.


Afterwards, the conditioned aerosol passes the measuring cell of the aerosol spectrometer.

The measurement zone is illuminated by light of two different wavelengths, which originates from two laser diodes with high lifetime. The light, which is scattered by the particles or droplets, is collected and recorded by a photodetector.



## Specifications

The optical signals are converted into electrical ones and are processed. Signal evaluation is performed by the firmware. Results like the particle size distribution and the particle number concentration are visualised on a display.

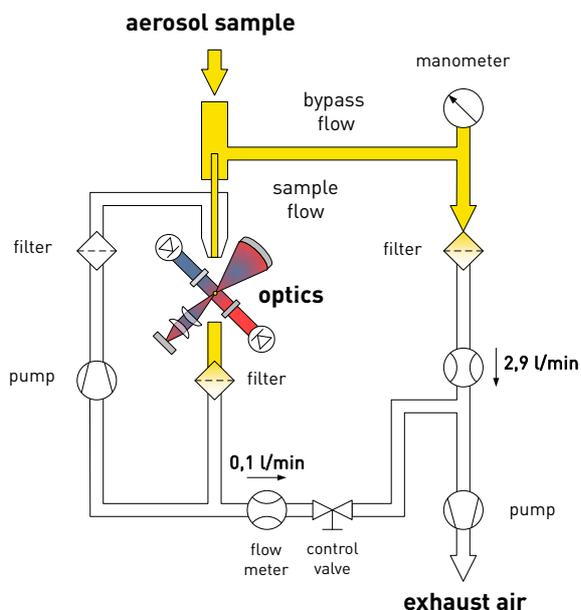

Principle of operation of the laser aerosol particle size spectrometer LAP 325.

After the analysis, the excess flow rate is continuously re-fed into the test channel. The excess flow re-feeding guarantees a constant flow rate through the test channel. Accordingly, aerosol sampling does not affect the pressure drop over the test specimen.

The implemented dilution is based on a closed, static dilution principle. For the FMS 375, the principle was automated and optimized regarding pressure-resistance. No additional dilution air is required.

By weighing out the good accessible absolute filters of the bypass flow and the sample flow path of the aerosol spectrometer, an additional gravimetric evaluation of the aerosol sample is possible.

The FMS 375 can be operated as stand-alone device via the firmware and the integrated display.

The instrument can also be operated by a computer and the software FMS375Win.

### Application example

The FMS375 is typically used for the characterisation of fractional separation efficiency of separators in engine or oil mist separation test systems.

During testing, an oil-containing droplet aerosol is fed into a flow channel. The test aerosol is analysed down and upstream of the oil mist separator.

The fractional separation efficiency is computed from the up- and downstream particle concentration ratio.

### Accessories

- pneumatic hose with quick coupling adaptor (DN 5)
- trolley kit (optional)

### Technical specifications

sample flow rate	3 l/min
analytical flow rate	0,1 l/min
size range	0,2 ... 40 µm
concentration range	< 4 10 <sup>8</sup> cm <sup>-3</sup>
aerosol inlet	8 mm (outer diameter)
temperature range	max. 120 °C
pressure range (inlet)	max. ± 800 mbar
power supply	110...240 V, 50...60 Hz
compressed air supply	min. 4 bar max. 8 bar
data interface	ethernet, USB
dimensions (w × h × d)	
w/o grip	530 × 440 × 500 mm
with grip	650 × 570 × 500 mm
weight (w/o accessories)	33 kg
normative references	ISO 17536 series of standards; VDI 3867-4:2020 (E); ISO 21501-1:2009

© Copyright 2020 Topas GmbH. Specifications are subject to change without notice.



QMS certified according to DIN EN ISO 9001.



12 100 11908 TMS

Topas GmbH  
Technologie-orientierte  
Partikel-, Analysen- und Sensortechnik  
Gasanstaltstraße 47 · D-01237 Dresden

Telefon +49 (351) 21 66 43 - 0  
Fax +49 (351) 21 66 43 55  
E-Mail office@topas-gmbh.de  
Internet www.topas-gmbh.de

TOPAS-GMBH DE

PARTICLE UNDER CONTROL